


Quick Manual JETI LiVal

Connect your *JETI* device to a PC or Laptop via USB or Bluetooth.

Start *JETI LiVal* under Programs in the Start menu (software and driver installation is described in the operation manual).

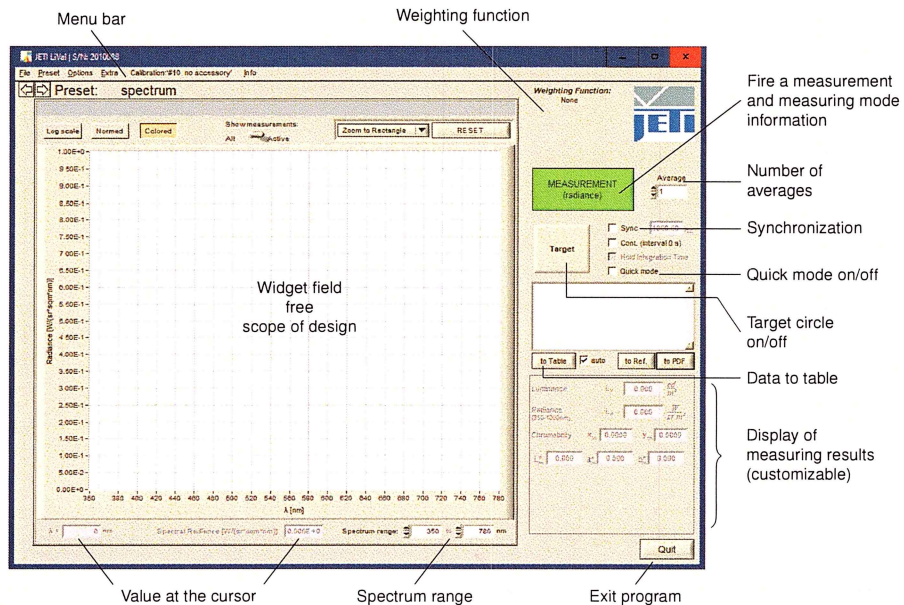

Measuring Mode Luminance/ Radiance

Attention: Do not look into the laser directly or with optical devices as telescopes or collimator lenses! Do not direct it to another person!

If *JETI LiVal* was started with no accessory attached to the device the software will be in Radiance mode.


Click the **Target** button and adjust the instrument viewing to the desired position.

Click the **Measurement** button to start the measurement. Maximum measuring time = 120 s.

Press **Esc** to stop a measurement in case of problems.

Spectrum and measuring values are shown.


If you need other measuring values, use the **Customize** option (right mouse click in measuring results field) or the **Preset** menu.

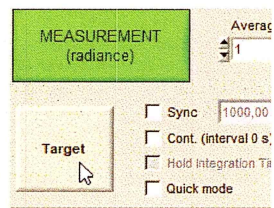

Quick Manual JETI LiVal

Connect your *JETI* device to a PC or Laptop via USB or Bluetooth.

Start *JETI LiVal* under Programs in the Start menu (software and driver installation is described in the operation manual).

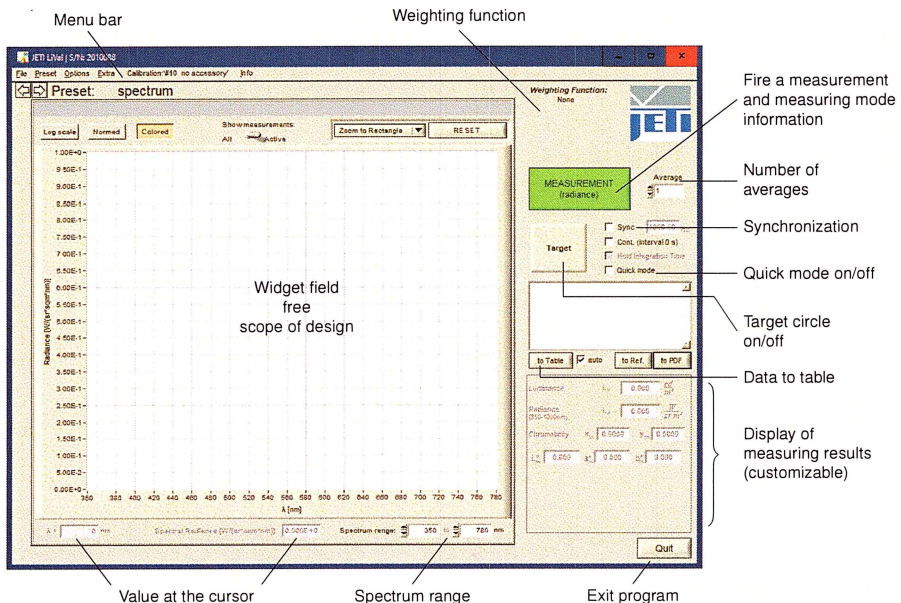

Measuring Mode Luminance/ Radiance

Attention: Do not look into the laser directly or with optical devices as telescopes or collimator lenses! Do not direct it to another person!

If *JETI LiVal* was started with no accessory attached to the device the software will be in Radiance mode.

Click the **Target** button and adjust the instrument viewing to the desired position.

Click the **Measurement** button to start the measurement. Maximum measuring time = 120 s.

Press **Esc** to stop a measurement in case of problems.

Spectrum and measuring values are shown.

If you need other measuring values, use the **Customize** option (right mouse click in measuring results field) or the **Preset** menu.

